



## PRODUCT STATEMENT

Product: PURUS™ Cleanroom Paper

### Autoclave Recommendations

PURUS cleanroom paper may be autoclaved with standard wet process provided the following procedures are taken:

1. The paper has been removed from its original cleanroom packaging.
2. The paper is in stacks no greater than 20 sheets for 22lb (80 gm) or 12 sheets for 28 lb (105gm).
3. Autoclave should not exceed 30 minutes of exposure at 121° C. Any variation of time, temperature or standard may be acceptable but is at the sole discretion of the user.
4. Paper should not be printed on prior to autoclave. If paper has been printed on, the ink used for printing should be capable of withstanding high temperatures. Wax based ribbons, inks or toners are NOT recommended. A high percentage resin-based ink is preferred.