

Non-Sterile Cleanroom Gowning Procedure

Step 1: Pre-entry



Bouffant/Hair Net Be sure to contain all hair.



Shoe Covers Contain all laces and tassels.

В



STEP 2: GOWNING



Gowning Gloves Recommended for Class 100 and better.



В



Mask Worn over or under hood. Bend nosepiece first for a snug facial fit.





Hood Ensure snug fit and proper face/neck seal.



STEP 3

Coverall

Step into the coverall. Be sure sleeves and upper garment do not

touch bench or floor.

Tuck shoulder panels from hood inside coverall before zipping up.




STEP 4

Boot Covers

Put on boots and pull boots (high-top shoe covers) over legs of coverall.

If cross-over bench is used, transfer each foot to cleaner side of bench as boot is donned.

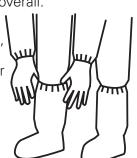

STEP 5

Goggles

Wear goggles/safety glasses when eye protection or additional particulate control is desired.



Step 6

Cleanroom Gloves (optional)

Place hem of glove over cuff of sleeve. Put on over gowning gloves or remove gowning gloves and then put on cleanroom gloves.



Non-Sterile Cleanroom Gowning Removal Procedure

STEP 1: IN ORDER



Gloves

If two pairs of gloves are worn, remove top pair of gloves, then discard.









Goggles/Safety Glasses Remove and properly contain.





Boot Covers Usually discard




after each change.



STEP 2

Coverall

If coverall is to be used again, either hang it in a controlled environment or prepare for storage.

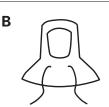

Hood

Hang in controlled environment.






C

Mask

Remove and usually discard.




STEP 3

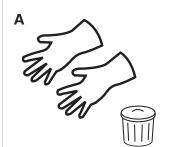


Gloves

It is recommended that gloves be discarded after each change.

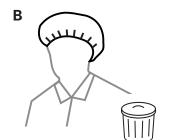

Bouffant/Hair Net Usually discard after each change.

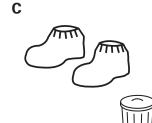

Shoe Covers Usually discard after each change.



CONTACT CLEANROOM CONNECTION FOR ALL OF YOUR CLEANROOM APPAREL. OUR **CUSTOMER SERVICE TEAM CAN HELP YOU** CHOOSE ALL OF YOUR CLEANROOM APPAREL.

CALL US AT 800.616.-5319

VISIT US ONLINE CLEANROOM SUPPLIES.COM



Sterile Cleanroom Gowning Procedure

STEP 1: PRE-ENTRY



Bouffant/Hair Net Be sure to contain all hair.




Shoe Covers Contain all laces

and tassels.

В



STEP 2: GOWNING

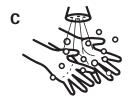

Wash hands using bactericidal soap and dry thoroughly.



First Sterile Gloves

Follow aseptic technique by touching only the inside of gloves. Sanitize.






Sanitize gloves after donning each article, if required.



STEP 3

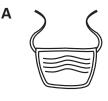

Mask

Adjust for a snug facial fit. May be worn over or under proper neck seal. hood.



В

Hood

Ensure snug fit and



STEP 4

Coverall

Carefully remove coverall from package by holding the inside neck area.

Unfold garment while preventing it from touching the floor, your clothes or bench.

Please exercise care when inserting leg to ensure donning snaps (if used) release and foot clears elastic ankle prior to transferring weight.



Don using aseptic technique.





STEP 5


Boot Covers

Put boots (high-top shoe covers) on using aseptic technique. Tuck coverall legs into boots.

The first boot donned can now be placed in the "clean" side of gowning area.



Don second boot and place the other foot in the "clean" side of the gowning area.

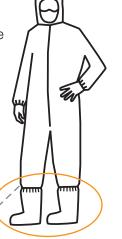

STEP 6

Goggles

Adjust for a snug facial fit.

Second Sterile Gloves Complete gowning steps

by aseptically donning second pair of sterile gloves over the first pair, making sure the cuffs are securely over the sleeves.



Perform a final inspection in gowning room mirror prior to entry into the sterile cleanroom.



Sterile Cleanroom Gowning Removal Procedure

REMOVAL OF STERILE GARMENTS

Enter gowning area from cleanroom and head directly to the "exit gown room" (preferred method).

If no "exit gown room" is available, go directly to the "non-clean" side, usually divided by a bench and/or line. Remove items in reverse order of gowning procedure and discard.








C



D

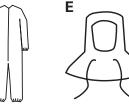







Outer Gloves

Goggles

Boot Covers

Coveralls

Hood

Mask

Inner Gloves Bouffant/Hair Net Shoe Covers

